

CytoflexSRT Filter Setup

Filter Position	Bandpass	Fluorochromes
VIOLET LASER (405nm)		
VioletSSC	405/10	
450Violet	450/45	CFP, e450, BFP, BV421, DAPI
525Violet	525/40	L/D AQUA, BV510
610Violet	610/20	BV605
660Violet	660/19	
780Violet	780/60	BV786 (BV785)
YELLOW LASER (561 nm)		
610Yellow	610/20	mKate, mRuby, PE-TexRd, mCherry
585Yellow	585/42	PI, PE, DsRed, mTomato, Miami Yellow
690Yellow	690/50	PE-Cy5
710Yellow	710/50	
780Yellow	780/60	PE-Cy7
RED LASER (638 nm)		
660Red	660/20	APC, Alexa Fluor 647, Cy5, DRAQ5
712Red	710/25	Alexa Fluor 700
780Red	780/60	APC-Cy7, APC-e780, APC-H7
BLUE LASER (488 nm)		
SSC	488/8	SSC
525Blue	525/40	FITC, eGFP, Alexa Fluor 488, YFP
690Blue	690/50	7AAD, PerCP-Cy5.5